CertoClav Connect

Laborautoklav

(User manual English: page 39)

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Laboratory autoclave

19 Table of contents

19	Table of contents.....	40
20	Preliminary note	42
21	CE - Declaration of Conformity	43
22	General safety instructions.....	44
23	Substances that must not be autoclaved.....	45
23.1	<i>Explosive substances.....</i>	45
23.2	<i>Flammable substances</i>	45
23.3	<i>Oxidizing agent.....</i>	45
23.4	<i>Combustible substances</i>	46
24	Autoclave - safety devices	46
25	Technical data	47
25.1	<i>Heating times.....</i>	47
25.2	<i>Cooling / evaporation times.....</i>	47
25.3	<i>Correlation table</i>	47
25.4	<i>CertoTouch technical data</i>	48
25.5	<i>Signs</i>	48
26	Scope of delivery.....	49
26.1	<i>The scope of delivery includes:</i>	49
26.2	<i>In addition, we offer the following accessories:.....</i>	49
27	Sterilization programs.....	49
27.1	<i>Standard programs for autoclaves with media sensor.....</i>	49
27.2	<i>Standard programs for autoclaves without media sensor.....</i>	50
28	Lineup	50
28.1	<i>Location.....</i>	50
28.2	<i>Autoclave.....</i>	50
28.3	<i>CertoTouch operating unit.....</i>	51
29	Operation	51
29.1	<i>Creating a user account on the CertoTouch</i>	51
29.2	<i>Before 1st use.....</i>	53
29.3	<i>Loading the device.....</i>	53
29.4	<i>Close lid.....</i>	54
29.5	<i>Start program.....</i>	55
29.6	<i>Venting</i>	55
29.7	<i>Sterilize / Autoclave</i>	55
29.8	<i>Status display</i>	56
29.9	<i>End of the cycle - cooling</i>	56
29.10	<i>Open and unload device</i>	56
29.11	<i>Cycle abort and power interruption.....</i>	57
30	Flexible media temperature sensor	57
30.1	<i>General.....</i>	57
30.2	<i>Application.....</i>	57
30.3	<i>Activating/deactivating the media sensor.....</i>	57
31	Maintenance and care of the autoclave.....	58
31.1	<i>Condensate tank.....</i>	58
31.2	<i>Cleaning of the device.....</i>	58
31.3	<i>Sealing ring.....</i>	58
32	Error messages	58
33	CertoTouch - control unit.....	59
33.1	<i>Support menu.....</i>	59
33.1.1	<i>Tutorial videos.....</i>	60
33.1.2	<i>Send email to CertoSupport</i>	60
33.1.3	<i>TeamViewer remote support and screen sharing.....</i>	61
33.2	<i>Logging menu.....</i>	62
33.2.1	<i>View logs</i>	62
33.2.2	<i>Print protocols</i>	62
33.2.3	<i>Example expression</i>	63
33.3	<i>Settings</i>	64

33.4	<i>User management</i>	64
33.5	<i>Network settings</i>	64
33.5.1	Connection to a WLAN network.....	64
33.5.2	Connection to a LAN network.....	65
33.5.3	Network communication	65
33.5.4	Connection of the autoclave to the local CertoServer.....	66
33.6	<i>Sterilization settings</i>	67
33.6.1	Export of all logs to USB stick.....	67
33.6.2	Export of all logs to SD card	67
33.6.3	Export of all protocols via network	68
33.6.4	Label Designer	68
33.7	<i>System settings</i>	69
33.7.1	Install updates via Internet.....	69
33.7.2	Install updates via USB stick	69
33.7.3	Install updates via SD card.....	70
33.7.4	Restore factory settings.....	70
33.7.5	Change language.....	71
33.7.6	Configuration of SMS and e-mail notification.....	71
33.8	<i>Sensor calibration</i>	72
33.8.1	Audit Trail.....	72
33.8.2	Lock settings	72
34	PC software	73
34.1	<i>CertoCloud Software Features</i>	73
34.2	<i>Display of sterilization protocols via web browser</i>	75
35	Transport and storage	76
35.1	<i>Storage conditions</i>	76
35.2	<i>Transport conditions</i>	76
36	Warranty	76

20 Preliminary note

Congratulations on your purchase of the CertoClav Connect small steam sterilizer.

CertoClav Connect is the most compact, fully automated laboratory autoclave on the market. It is operated by means of a 7 inch touch screen. Operating steps are explained with videos directly on the touch screen. The software includes user management, is GLP/GMP compliant and allows paperless documentation. Individual programs can be created directly on the screen.

If the device is connected to the Internet via WLAN or LAN, the autoclave is connected to the CertoCloud via a secure connection. This allows all protocols to be accessed browser-based, and individual programs to be created and sent to devices. Programs created on the device are also synchronized.

Important features

- Fully automatic autoclave
- Suitable for sterilization of waste, culture media, liquids and simple solids
- 18 liters chamber volume
- Temperature programmable from 25°C up to a maximum of 140°C
- Maximum 14 days sterilization time at a time
- Automatic venting
- Electric exhaust valve
- Optional flexible media sensor
- Time and temperature can be selected individually
- Controller with LCD display
- Venting in the fractional flow process
- Ethernet, LAN, WiFi
- 7" tablet with video assistant
- Cloud connection
- Heating integrated in the floor
- Electronic safety interlock

Contact by e-mail: support@certoclav.com

Your CertoClav Team

21 CE - Declaration of Conformity

This is to certify that the **CERTOCLAV CONNECT** autoclave complies with the following EU directives:

2014/35/EU, Low Voltage Directive

Related standards:

- EN 61010-2-040:2015 Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 2-040: Particular requirements for sterilizers and washer-disinfectors for the treatment of medical material
- EN 61010-1:2010 Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements

2014/35/EU, Low Voltage Directive

2014/30/EU, EMC Directive.

Related standards:

- EN 61326-1:2013 Electrical equipment for measurement, control, and laboratory use - EMC requirements - Part 1: General requirements (intended date of replacement: EDIN EN 61326-1:2018)
- EN 61000-3-2:2014 Electromagnetic compatibility (EMC) - Part 3- 2: Limits - Limits for harmonic current (equipment input current ≤ 16 A per phase)
- EN 61000-3-3:2013 Electromagnetic compatibility (EMC) - Part 3- 3: Limits - Limitation of voltage changes, voltage fluctuations and flicker in public low-voltage distribution systems for equipment with a rated current ≤ 16 A per conductor and not subject to special connection conditions

2011/65/EU Directive on the restriction of the use of certain hazardous substances in electrical and electronic equipment

Related standards:

- EN 50581:2012 Technical documentation for the evaluation of electrical and electronic equipment with regard to the restriction of hazardous substances.

Manufacturing quality:

CertoClav manufactures according to modern manufacturing methods using high quality materials. Decades of experience and proven employees ensure high quality products and maximum customer service. All devices are subject to 100% testing for function and quality. CertoClav products are manufactured exclusively in Austria.

Manufacturer:

CERTOCLAV Sterilizer GmbH
A-4060 Leonding
Austria

Michael Dirix
Managing Director

Leonding, January 2021

22 General safety instructions

- ⚠ Use the autoclave and accessories only in accordance with the operating instructions
- ⚠ Keep the operating instructions in a place that is accessible to all users of the device.
- ⚠ Only use the supplied screen to control the autoclave.
- ⚠ Never tamper with safety equipment. Always use the autoclave and control of the same series. Do not use lids and chambers of different series together.
- ⚠ The surface of the device (metal and plastic) becomes hot during operation. Suitable protective gloves must be used.
- ⚠ This device must not be used in potentially explosive atmospheres.
- ⚠ Do not mount the device near high power electronics.
- ⚠ The CertoTouch screen must not be closer than 30cm to a CertoClav autoclave (heat protection).
- ⚠ Disconnect the power plug before opening the housing.
- ⚠ Only use the device with the intended mains voltage.
- ⚠ Do not load, bend or twist the cables. Do not place heavy objects on the cables.
- ⚠ Intended for laboratory use by trained personnel only. Do not use in the household. Not suitable for medical purposes.
- ⚠ Never operate the device without water.
- ⚠ Target level: 1.5 liters (minimum 0.7 liters, maximum 2.5 liters).
- ⚠ Do not sterilize pressure-tight sealed vessels or packages, as there is no pressure equalization.
- ⚠ Use only water to generate steam.
- ⚠ Only fill bottles to a maximum of 75% with liquid to prevent boiling over.
- ⚠ Do not release steam in the direction of the person - risk of burns. For this purpose, use the supplied steam exhaust container on the backpack of the unit!
- ⚠ Wait until the program has ended before opening the device. If liquids are being sterilized, the autoclave can only be opened from 80°C media temperature.
- ⚠ When autoclaving liquids, there is a risk of boiling delay. To avoid this, the autoclave should only be opened if the media temperature is below 80°C.
- ⚠ Do not transport the autoclave in a filled state.
- ⚠ The device may only be repaired by CertoClav Sterilizer GmbH or by persons authorized by CertoClav Sterilizer GmbH (specialist company or customer service) using original spare parts and observing the service instructions.
- ⚠ Remove the power plug before unscrewing the electronics housing or the base.
- ⚠ Do not use force to open the lid under pressure. It is electronically locked.
- ⚠ Never sterilize flammable or explosive materials in the device.

- ⚠ In case of damage to the electrical supply line, the device must be disconnected from the mains immediately.
- ⚠ Never insert objects into the ventilation slots. Do not pour any liquids into the ventilation slots. Damage due to electric shocks and short circuits.
- ⚠ To avoid corrosion, do not sterilize substances with a high salt content.
- ⚠ Do not empty the condensation tank until it has cooled down - risk of scalding

23 Substances that must not be autoclaved

Never use the autoclave to sterilize any of the following materials or alkaline solutions. Sterilization of such substances may cause an explosion, corrosion of the sterilization chamber, piping and wear of the seals.

23.1 Explosive substances

- Nitroglycol, nitroglycerin, nitrocellulose and other explosive nitric acid esters.
- Trinitrobenzene, trinitrotoluene, picric acid and other explosive nitro compounds.
- Peracetic acid, methyl ethyl ketone peroxide, benzoyl peroxide and other organic peroxides.

23.2 Flammable substances

- Metallic lithium, potassium, sodium, yellow phosphorus, phosphorus sulfide and red phosphorus.
- Celluloids, calcium carbide (carbide), calcium phosphide and magnesium powder.
- Aluminum powder, magnesium powder and metal powder except aluminum powder.
- Sodium dithionite (or sodium hydrosulfite).

23.3 Oxidizing agent

- Potassium chlorate, sodium chlorate, ammonium chlorate and other chlorates.
- Potassium perchlorate, sodium perchlorate, ammonium perchlorate and other perchlorates.
- Potassium peroxide, sodium peroxide, barium peroxide and other inorganic peroxides.
- Potassium nitrate, sodium nitrate, ammonium nitrate and other nitrates.
- Sodium chlorite and other chlorites.
- Calcium hypochlorite and other hypochlorites.

23.4 Combustible substances

- Ethyl ether, gasoline, acetaldehyde, propylene oxide, carbon disulfide and other substances whose flash points range from -30 to 0 °C.
- Methanol, ethanol, xylene, benzyl acetate (or amyl acetate) and other substances whose flash points are between 0 and 30 °C.
- Kerosene, gas oil, turpentine oil, isopentyl alcohol (or isoamyl alcohol), acetic acid and other substances with a flash point between 30 and 65 °C.
- Combustible gas (hydrogen, acetylene, ethylene, methane, ethane, propane, butane and other substances operating at a temperature of 60 °C at an absolute pressure of 0.2 bar)

If a saline solution has been spilled inside the chamber, remove all liquid from the chamber and wipe up the liquid around the lid seal. Clean the lid seal with distilled or demineralized water.

24 Autoclave - safety devices

Locking mechanism on the handle

The CertoClav Secure Lock locking mechanism secures the lid of the autoclave against opening under pressure. Turning the red lever to "Open" automatically opens the exhaust valve and releases steam. As long as the red lever is not set to "Closed", the exhaust valve remains open and prevents pressure build-up.

Locking mechanism on the lid

As soon as a program has been started or the temperature in the chamber exceeds 90°C or the temperature in bottles exceeds 80°C, the lid locks electronically.

Safety valve in handle

Both handles are equipped with a pressure relief safety valve

Safety valve on the lid

The safety valve on the lid opens from a pressure of 3.6 to 4 bar and closes again when the operating pressure is reached. In this case, replace the pressure relief valve, as there is probably a defect.

CAUTION: In this case, be sure to send the device in for inspection.

25 Technical data

Dimensions and weight			
Material	Stainless steel 304L	Certificates	CE
Dimensions			
Chamber volume	18 liters	Device height	60 cm
Chamber diameter	26 cm	Device width	44 cm
Chamber height	35 cm	Device depth	40 cm
Chamber usable height	30 cm	Device weight	15 kg
Operating data			
Max. Operating temperature	140 °C	Power supply	230V, 50/60 Hz
Max. Operating pressure	2.7 Bar (39 psi)	Power	1300 W
Test print	4.1 Bar (59 psi)		

25.1 Heating times

Loading quantity	Heating time from 25 °C to 121 °C
Low loading	15 - 20 minutes
High load	25 - 40 minutes

25.2 Cooling / evaporation times

Loading quantity	Cooling time at fast evaporation rate from 121 °C to 80 °C	Cooling time at slow evaporation rate from 121 °C to 80 °C
Low	20 - 35 minutes	40 - 60 minutes
High	40 - 60 minutes	80 - 120 minutes

25.3 Correlation table

T in °C	p in bar (kgf/cm ²)	T in °C	p in bar (kgf/cm ²)
100	0,03	118	0,90
101	0,07	119	0,96
102	0,11	120	1,02
103	0,15	121	1,09
104	0,19	122	1,15
105	0,23	123	1,22
106	0,28	124	1,29
107	0,32	125	1,37
108	0,37	126	1,44
109	0,41	127	1,52
110	0,46	128	1,59
111	0,51	129	1,67
112	0,56	130	1,75
113	0,61	131	1,84
114	0,67	132	1,92
115	0,72	133	2,01
116	0,78	134	2,10
117	0,84	135	2,19

25.4 CertoTouch technical data

Screen type	7" capacitive touch screen, LED backlighting
Screen resolution	800 x 480 pixel
Screen brightness	450cd/m ²
CPU	Freescale iMX.53 x 800MHz / 1.0GHz
Core	ARM Cortex-A8
Memory card connectors	Micro SD and USB stick optional
LAN	Ethernet 10/100M (RJ45) optional
W-LAN	IEEE 802.11
Accessories connections	CertoClav barcode scanner connector CertoClav label printer connection
Audio	Headphone jack Loudspeaker
Loudspeaker	1W
Power consumption	9W
Working temperature	-10°C to 65°C
Storage temperature	-25°C to 80°C
Power supply	DC 9V to 36V
Dimensions	220x132x36.5mm (length/width/height)
Weight	740g
Screen protection level	All-round splash water protection IP44

25.5 Signs

	Directive 2012/19/EU (WEEE): The device must not be disposed of with residual waste, but only by public waste disposal services and subsequent return to the manufacturer and importer.
	WARNING: Read the instructions before use
	WARNING: Hot surface
	The CE mark confirms the conformity of the device to the directives NSR 2014/35 EU + EMC 2014/30 EU, as well as EN 61010-1:2010.

26 Scope of delivery

After you have opened the packaging, please check the product for transport damage. If you find any damage, please contact your dealer and do not operate the device for safety reasons. Please also check whether the full scope of delivery is included.

26.1 The scope of delivery includes:

- Autoclave Connect including lid
- CertoTouch control panel with stand, 4 mounting screws and 12V power supply unit
- Interface box
- RS-232 connection cable
- Autoclave cold appliance power cable
- Instrument plate with water level indicator
- The CertoCloud basic version is available at www.certocloud.com.
- Instruction manual
- Warranty certificate
- Media sensor (optional depending on the device version)

26.2 In addition, we offer the following accessories:

- Art. No. 1300149 Wire basket with handle, \varnothing 25 cm, stainless steel.
- Art. No. 1310186 Service set 2 with all seals
- Art. No. 8500482 Water distiller. It is recommended to use distilled water
- Art. No. 1300187 Protocol printer, with connector plug.
- Art. No. 8500496 Thermal paper for log printer.
- Art. No. 8500508 CertoServer
- Art. No. 8500212 CertoCloud Premium Account

For more information about our assortment please visit: www.certoclav.com

27 Sterilization programs

27.1 Standard programs for autoclaves with media sensor

Name	Sterilization temperature [°C]	Sterilization time [min]
Solids 121	121	20
Solids 134	134	10
Solid 140	140	5
Liquid 121	121	20
Liquid 134	134	20
Liquid 140	140	20

Note: The programs can be changed as desired. See the chapter on editing programs.

27.2 Standard programs for autoclaves without media sensor

Name	Sterilization temperature [°C]	Sterilization time [min]
Solids 121	121	20
Solids 134	134	10
Solid 140	140	5
Liquid 121 1000ml	121	30
Liquid 121 3000ml	121	40
Liquid 121 5000ml	121	50
Liquid 140 1000ml	140	30
Liquid 140 3000ml	140	40
Liquid 140 5000ml	140	50

Note: The programs can be changed as desired. See the chapter on editing programs.

28 Lineup

28.1 Location


-  Avoid ambient temperatures below 5° and above 35 °C and humidity above 80 %.
-  Do not block the side ventilation slots of the device.
-  Avoid locations subject to shock and vibration and do not use the device outdoors.
-  Make sure that your instrument is not set up in the vicinity of steam-sensitive laboratory equipment (e.g. microscopes).

28.2 Autoclave

- Take the unit out of the box by the two handles and place it on a stable surface so that an ideal working height is provided.
- Remove all packing materials.
- Plug the power cord into the connector on the backpack of the device and connect it to a suitable power outlet with identical parameters as indicated on the nameplate of the autoclave.
- Switch on the device with the green rocker switch.
- The device is now automatically unlocked.
- Press down the red rotary knob on the right lid handle and turn it clockwise.
- Then turn the lid counterclockwise to open the autoclave.
- Remove the CertoTouch and the interface box from the chamber.
- If necessary, remove the accessories supplied.

28.3 CertoTouch operating unit


- Remove the device from the packaging.
- Use the 4 screws provided to secure the screen to the stand.
- Place the control unit next to the device. The stand can be used to set a suitable viewing angle.
- Connect the touch screen to the interface box using the cable provided.
- Connect the interface box with the power adapter to the power outlet.
- Connect the autoclave to the interface box by plugging the RS232 cable into the connector on the backpack of the device and into the COM4 position of the interface box.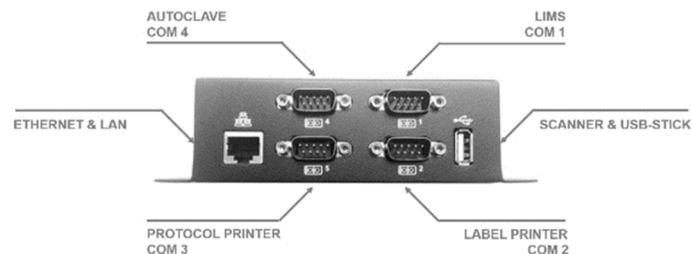


29 Operation

29.1 Creating a user account on the CertoTouch

To switch on the CertoTouch, press the -switch on the front of the CertoTouch for at least 3 seconds.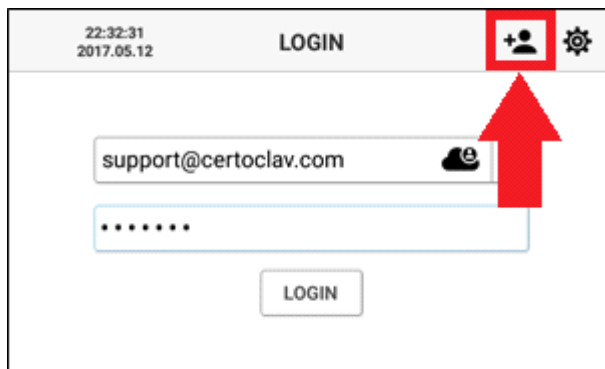

After switching on the touch screen, the login window appears. To create a new user, press the plus symbol at the top right edge of the screen.

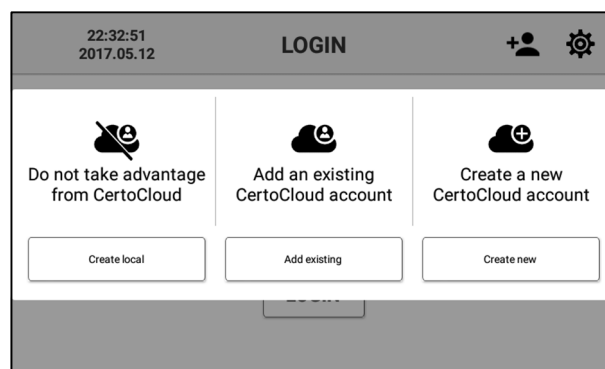To create a local account without server connection, press the "Local" button on the left side of the selection window. To add an existing CertoCloud user account with server connection press the middle button. If you want to create a new CertoCloud account with server connection, press the Create new button on the right side of the selection window.

To create an account, enter your user data. Text fields can be selected and edited by finger pressure. Enter your telephone number including country code. This is required if you want to receive status notifications of the autoclave via SMS. Enter a password of your choice with at least 4 characters repeatedly. To cancel the registration, press the Back button in the upper left corner of the screen. To complete the registration, press the Register button.

To add an existing CertoCloud account, enter your email address and password. Then press Add account. To cancel the registration, press the Back button in the upper left corner of the screen.

To create a CertoCloud account, enter your user data in the text fields provided. Please enter your cell phone number including country code. To confirm your email address and phone number, you will receive an email and SMS. Enter the password of your choice twice and then press the Register button.

To log in to a recently created account, first select your email address from the list and then enter your password. If you want to use a server-connected account, make sure that the online mode is enabled. To enable online mode, navigate to Settings->Device->Online Mode.

Activate max.wagner@mylab.com

Activation key

Activate account

Send activation key to max.wagner@mylab.com

If you are logging into your CertoCloud account for the first time, you will be asked for an activation key. You received the activation key via SMS and email after you created your CertoCloud account. Enter this key in the text field and press Activate account. Now your CertoCloud account is activated and you can now log in with your CertoCloud account.

29.2 Before 1st use

-  Read the operating instructions completely before using the autoclave for the first time.
-  Perform the first sterilization cycle without loading (with water only).

The first test run serves to check the function. It is intended to detect any damage that may have occurred during transport.

29.3 Loading the device

- Place the instrument plate in the chamber.
 - Fill the chamber with distilled water up to the indicator on the bottom plate. (min. 0.7, max. 2.5 liters)
 - The device can also be filled with a larger amount of liquid. However, this leads to long heating times and very slow cooling.
 - Place the load in the chamber. It is important that the steam can circulate freely, therefore the load must have a distance to the chamber wall!
 - It is recommended to use a wire basket for destruction bags or similar to avoid blocking the connections and to ensure sufficient steam supply.
 - Be careful not to overfill the chamber, otherwise the steam cannot circulate properly and the sterilization may be defective as a result!
 - When autoclaving liquids, do not fill containers fuller than $\frac{3}{4}$ of capacity to prevent boiling over.
 - When liquefying culture media, the vessel should not be filled with more than 2 liters.
-  **Do not autoclave tightly sealed vessels, as there is no pressure equalization.**


29.4 Close lid

- Check that the white lid gasket is properly inserted in the lid.
- Before closing, check that there are no foreign bodies between the underside of the lid and the seal, otherwise steam could escape.
- Be careful not to pinch your fingers when closing.


(Symbol image)

- Press the red rotary knob and turn it clockwise to the OPEN position.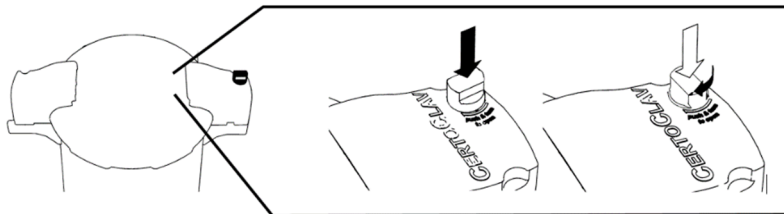


- Place the lid on the sterilization chamber and turn it clockwise until the handles are on top of each other.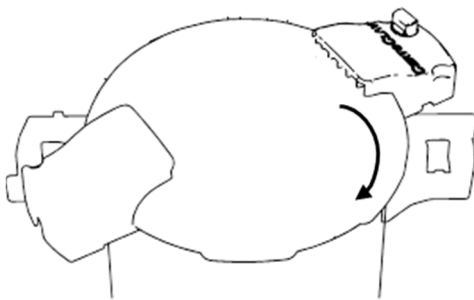

- Turn the red rotary knob counterclockwise to lock the lid.

29.5 Start program

- Check that the temperature in the chamber is below 50°C before starting the cycle.
- On the CertoTouch, select the 'Sterilization' tab to access the program overview.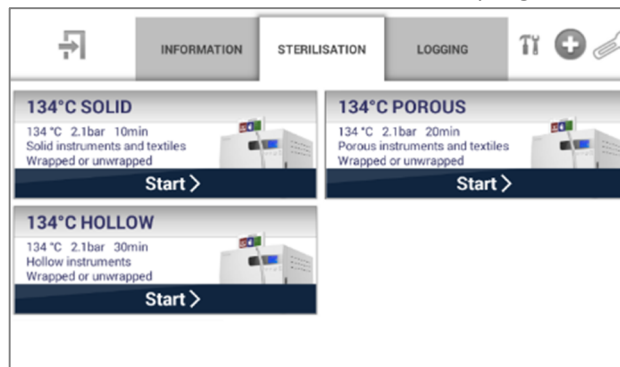

- Select the desired program and press Start.
- Follow the step-by-step videos if necessary. The program will start automatically afterwards.

29.6 Venting

If the "Venting" program option is activated, the chamber is vented at the beginning of the sterilization cycle at 100°C for 10 minutes. After venting, saturated water vapor remains in the chamber. The exhaust valve is then closed.

Venting of the autoclave is necessary to achieve uniform temperature distribution in the sterilization chamber and optimal temperature transfer.


29.7 Sterilize / Autoclave

The selected program is now executed. If the permissible temperature band of -1°C/+2°C deviates from the target temperature, the program is aborted!

- ⚠ When sterilizing Petri dishes or empty containers, the air remaining in the container expands and can possibly increase the pressure in the chamber extraordinarily. When the pressure exceeds the saturated steam pressure in the chamber, the automatic exhaust valve opens and allows the air to flow out of the chamber into the condensate container.

29.8 Status display

As soon as a program is running, the status screen is displayed. The status screen displays all relevant data about the current program.

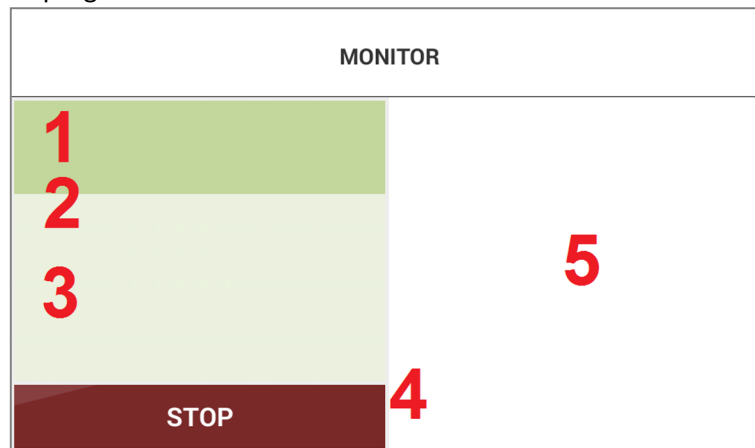- | | |
|----------|---|
| 1 | Name of the current program |
| 2 | Status of the running program. The status can be: "Program running", "Completed" "Not started", "Preparing to start". |
| 3 | A list of the program steps |
| 4 | Current temperature of sterilization chamber and media sensor in [°C].
Pressure in the sterilization chamber in [bar]. |
| 5 | Time counter and program history as graph. Alternatively as plain text. |

29.9 End of the cycle - cooling

- As the last step, the control unit cools down to 90°C or to 80°C in the liquids program. After this, a message appears on the display that the lid can be opened.

29.10 Open and unload device

In order to be able to open the device and remove the items to be sterilized, the chamber temperature must have dropped to 80°C. Only then does the electronic lock open and the lid can be opened. This process serves to protect the personnel from scalding when opening the lid.

- Turn the red knob to open, which opens the exhaust valve. Any residual steam escapes.
- Turn the lid counterclockwise, remove it and swing it slightly to allow residual water to drain.
- To improve drying, remove the basket from the chamber and allow it to cool to room temperature before removing the load.

- ⚠ When opening the lid after sterilization, never hold your face or hands over the open chamber. Risk of scalding due to escaping steam.
- ⚠ It is best to open the lid with the handle only a few centimeters at first so that the first steam can escape. To do this, stand at a sufficient distance from the device. Then carefully open the lid completely.
- ⚠ Always use heat-insulating protective gloves to remove materials from the chamber.
- ⚠ Due to the boiling delay, liquids need a lot of time to cool down. Make sure that the temperature of the liquid has dropped sufficiently before removing it from the chamber.
- ⚠ **INCOMPLETE CYCLE:** If, for whatever reason, the cycle is interrupted (by switching off or power failure), sterilization has not been performed properly and must be repeated. Before opening, follow the safety instructions for cooling.

29.11 Cycle abort and power interruption

- ⚠ If the cycle is interrupted during the sterilization of liquids, the lid will only be unlocked electronically once the chamber temperature is below 80 °C. The cycle must be restarted.
- ⚠ In the event of a power failure, the lid remains locked. For safety reasons, the device cannot be opened when it is switched off. To do this, the device must first be supplied with voltage again and switched on. Only then will the electronic locking mechanism open.

30 Flexible media temperature sensor

30.1 General

The flexible media temperature sensor is used to measure the temperature in the medium to be sterilized. Since boiling distortion occurs especially during the sterilization of liquids, it makes sense to use the flexible media sensor. In addition, after the sterilization cycle, you can check exactly which temperatures the medium was exposed to.

30.2 Application

After loading the sterilization chamber, place the media sensor in a reference container that corresponds to those of the sterilized material. If possible, place the reference jar in the center of the chamber and ensure that the sensor is as centered as possible in the medium.

30.3 Activating/deactivating the media sensor

You can define for each program in the program editing whether the medium sensor is to be activated for a specific program. With an activated media sensor, the temperature in the medium is controlled. Only when the medium has reached the set temperature does the holding phase begin. The sterilization time thus does not start until the medium has also reached the sterilization temperature.

- ⚠ Do not twist, knot or pull the cable of the temperature sensor. A damaged cable may cause malfunction or failure of the sensor.
- ⚠ Make sure that the temperature probe does not come into contact with soft components of the Petri dishes when sterilizing plastic Petri dishes. When the temperature has dropped, the plastic hardens and it may not be possible to remove these components from the temperature probe

31 Maintenance and care of the autoclave

31.1 Condensate tank

Water collects in the condensate container during each sterilization process. Please empty the condensate container after each cycle. Before each new cycle, please check whether the condensate container has been emptied.

- ⚠ Do not empty the condensate container before the autoclave has cooled down to 40°C.
- ⚠ If the condensate container has not been inserted, the exhaust valve is blocked.

31.2 Cleaning of the device

Do not use aggressive cleaning agents to clean the device housing and use a soft cloth to avoid scratching the housing and chamber.

31.3 Sealing ring

A sealing ring that has become brittle at the edge over the course of time could cause an outflow of the cause steam. The average service life is 2 years. Depending on the number of operating hours, this period may vary.

32 Error messages

Should a critical malfunction occur in the autoclave, for safety reasons a Error message triggered. The error message does not disappear until it is confirmed.


Error code	Malfunction	Measure
Error 1	Fan defective	Replace fan

Error 2	Lid is not closed	Please close the lid before starting a program. Follow the displayed step-by-step video for this.
Error 3	Water level in the chamber too high	A maximum of 2.5 liters of water should be in the chamber before a sterilization cycle. Check that the amount of water is correct.
Error 4	Water level in the chamber too low	Please fill at least 0.7 liters of water into the sterilization chamber before starting a program.
Error 5	Lid does not lock	Please turn off the device and check if something is blocking the mechanism. Then try it again. If the error message is still displayed, contact the service personnel.
Error 6	Timeout during heating process	Target temperature cannot be reached. The autoclave may be empty or the heater may be defective. Please contact the service staff.
Error 7	Maximum permissible pressure exceeded	Please contact the service staff.
Error 8	Maximum permissible temperature exceeded	Possible errors: Heating defective or temperature sensor defective. Please contact the service personnel.
Error 9	Pressure not in the permitted range	Unexpected pressure fluctuation in the chamber. Please do not place tightly closed vessels in the sterilization chamber.
Error 10	Chamber temperature not within permissible temperature range	Please contact the service staff. Temperature fluctuations during sterilization. Possible cause of error: heating defective, temperature sensor defective
Error 11	Media temperature not within the permissible temperature range	Please contact the service staff. Temperature fluctuations during sterilization. Possible cause of error: Temperature sensor defective
Error 13	Temperature band not maintained	Please contact the service personnel. Collective error of Error 10 and Error 11
Error 14	Cycle aborted due to a critical error	If a critical error occurred during sterilization and a sterilization process had to be aborted as a result, Error 14 is also displayed
Error 15	Cycle aborted due to user interaction	No action required. The program was canceled manually. Please wait until the chamber temperature has dropped below 80 °C before opening the unit
Error 16	Connection failure between tablet PC and autoclave	Please check the cable connections between the device, interface box and CertoTouch control unit.

33 CertoTouch - control unit


33.1 Support menu

Navigate to the "Information Tab" to display the information menu. Here it is possible to view training videos and contact CertoSupport directly from the touch screen.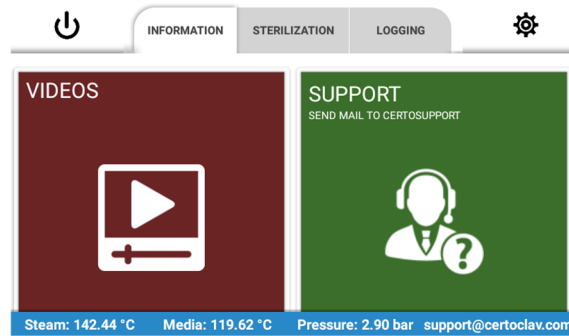


33.1.1 Tutorial videos

To scroll through the list of all available videos, press the Video section in the Information tab. Select the video you want to watch in full-screen mode.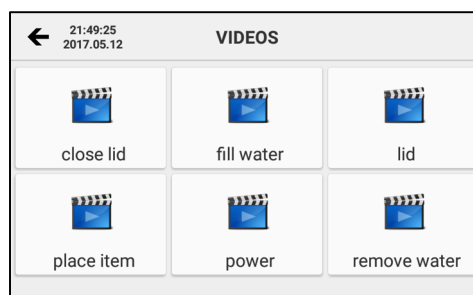

33.1.2 Send email to CertoSupport


To send an e-mail to CertoSupport directly from the device, the device must be connected to the Internet. To connect the device to the Internet, you can plug a network cable into the Ethernet port or establish a WLAN connection to your WLAN router. Press the "SUPPORT button" to open the email window.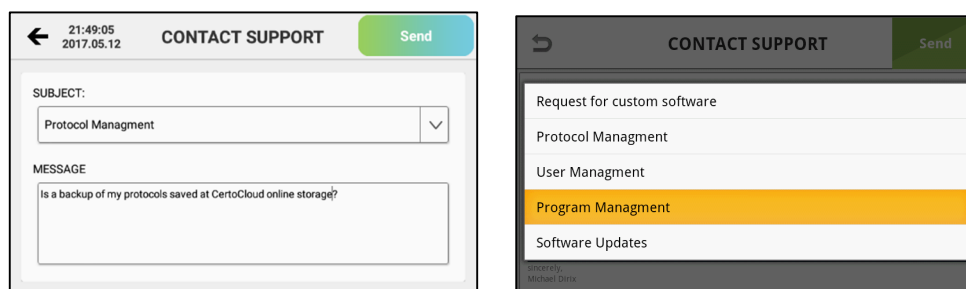


Select the subject of your message to CertoSupport from the drop-down list. Then enter your message to CertoSupport in the message text field. To send the e-mail, press the Send button.

33.1.3 TeamViewer remote support and screen sharing

To share the screen with CertoSupport using TeamViewer, navigate to the Information tab and press the Support button. Remote support is only available by prior appointment via email or phone. As soon as you start a TeamViewer session, a window will appear displaying your TeamViewer ID. Give this ID to the remote support team via phone or email.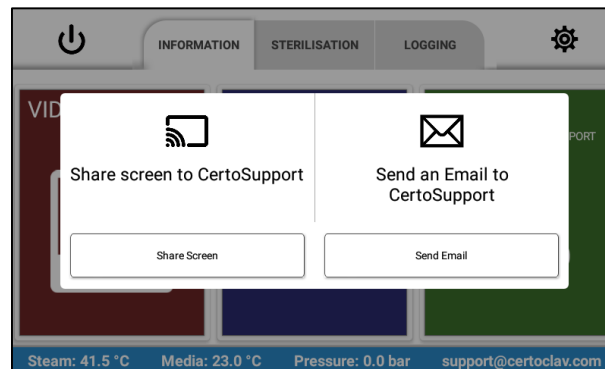


Wait until the support team joins the TeamViewer session. Confirm the incoming connection by pressing the Allow button. Now the technician can access the menu of your device. To cancel the screen transfer, press the X icon on the screen.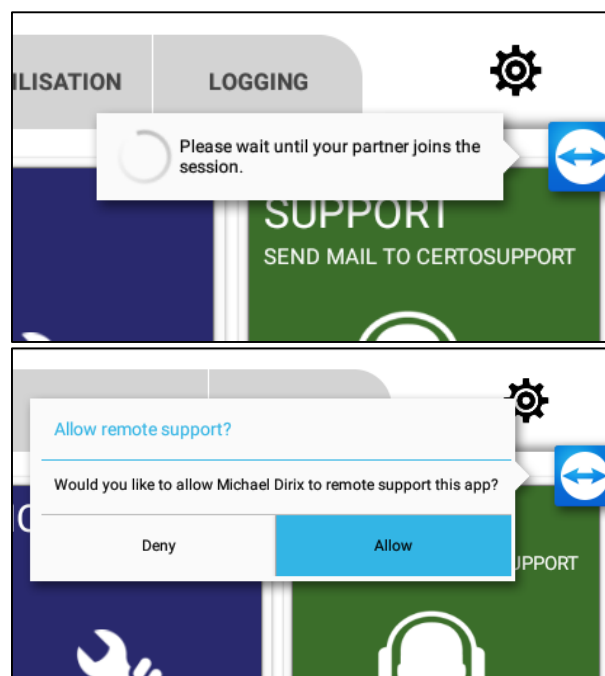

33.2 Logging menu


33.2.1 View logs

To view sterilization logs, navigate to the Logs tab. A list of all recorded logs is displayed on the left half of the screen. The cloud icon in the log list indicates whether the logs have been saved in the CertoCloud or on your local CertoServer. Press the drop-down menu above the log list to sort the logs by specific criteria. To view the details and graphical view of a log, select the log with a finger press on the list item. The graphical history of the log is now displayed on the right side of the window. To display the log as a list, press the list icon above the graph.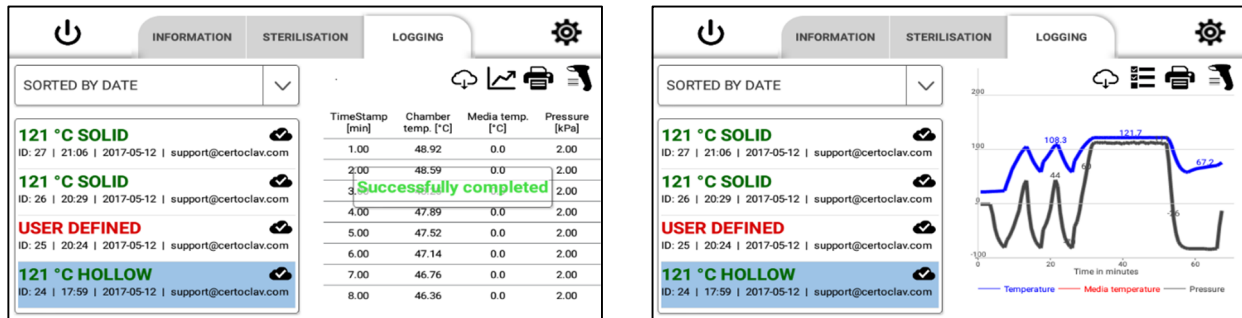

33.2.2 Print protocols

- 6) Make sure that a log printer is connected to the device.
- 7) Navigate to the Logging tab.
- 8) On the left side of the screen, select the protocol you want to print.
- 9) Press the print icon at the upper right corner of the screen.
- 10) Select the Print log option

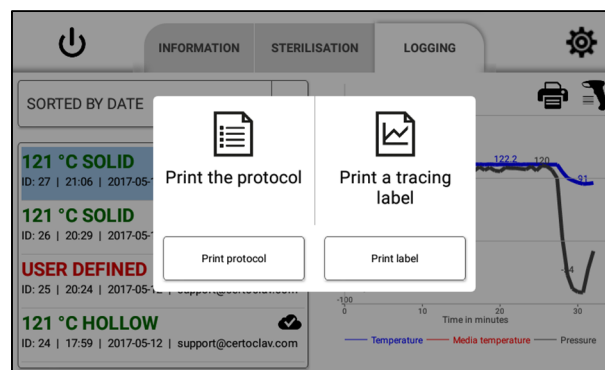33.2.3 Example expression

```
Header line 1
Header line 2
Header line 3
Model:      CertoClav Connect
S/N.:       C01.0001
Program:     121°C LIQUID
Program description:
20 minutes at 121°C
0 minutes at 60°C
Media:      NO
Fast Cooling:NO
Cycle no:    971
Start time: 12.01.15 17:42:00
End time:   12.01.15 18:16:00
Result:     PASSED
hh:mm:ss    Temp °C    Press [bar]
17:42:00    27.5           1.00
17:42:20    27.5           0.50
17:42:40    27.6           0.18
17:43:00    80.4           0.80
[...]
18:15:00    65.9           0.18
18:15:20    65.4           0.18
18:15:40    68.0           0.50
18:16:00    70.2           1.00
Program finished successfully
END

Signed by _____
Confirmed by _____
```

Optional

33.3 Settings

Press the  icon at the top right of the screen to enter the settings menu, where the following settings can be made:

33.4 User management


On the right side of the window all registered users are now displayed. Here you can change the user data and passwords with the Edit button and remove a user account from the touch PC with the Delete button.

33.5 Network settings


33.5.1 Connection to a WLAN network

4. Click on the Network entry.
5. If the WiFi checkbox is not set, press the empty checkbox to enable WiFi.
6. Then press Manage WLAN Connections to display a list of available WLAN networks. Select the relevant WLAN network and enter access keys and other parameters if necessary.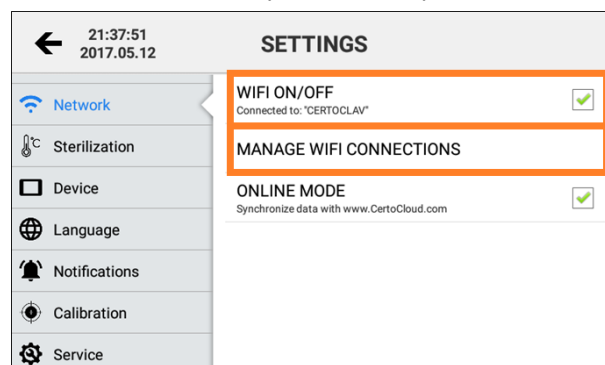

33.5.2 Connection to a LAN network


The autoclave can also be connected to a LAN network using an Ethernet cable. Connect the network cable to the Ethernet socket of the interface box.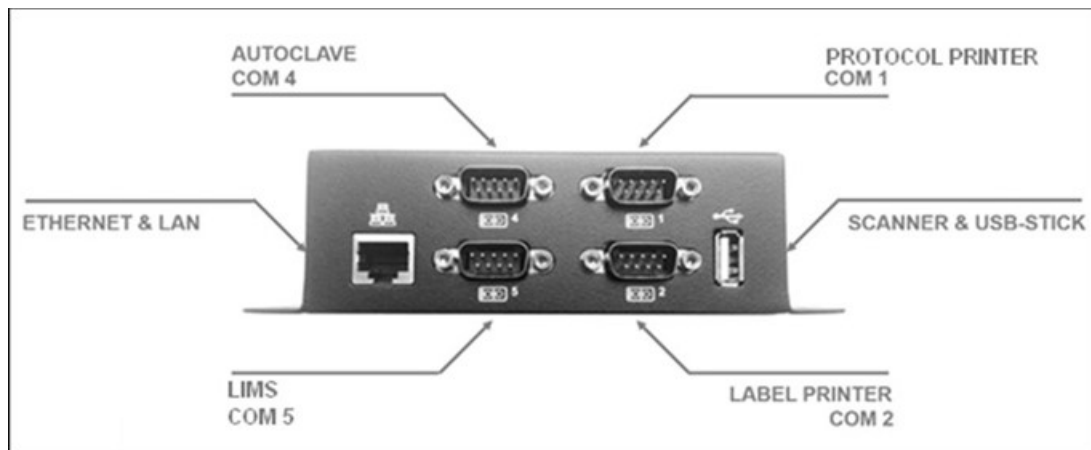

33.5.3 Network communication


If the data of the autoclave is to be synchronized with the CertoCloud (www.certocloud.com) or the local CertoServer (<http://certocloud>), the online mode must be activated.

3. Select the Network entry on the left side of the settings window.
4. Press the Online mode entry to activate or deactivate data synchronization with CertoCloud. Note: If the online mode is deactivated, your logs are only saved locally on the autoclave. As soon as you reactivate the online mode, all logs saved locally until then will be uploaded to CertoCloud and are thus also available online in your backup storage.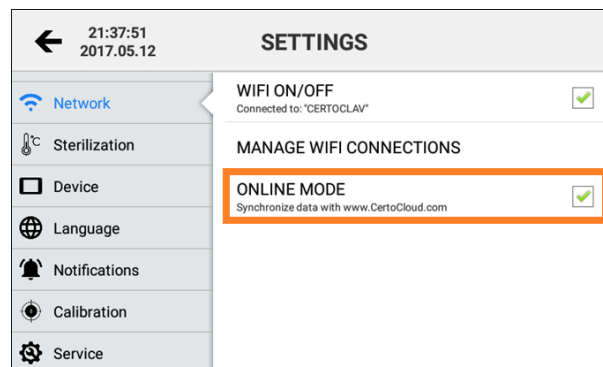


33.5.4 Connection of the autoclave to the local CertoServer


7. Connect the CertoServer to the interface box or the in-house network.
8. Connect the interface box to the same network as the CertoServer.
9. Navigate to the button: Settings->Network->Select server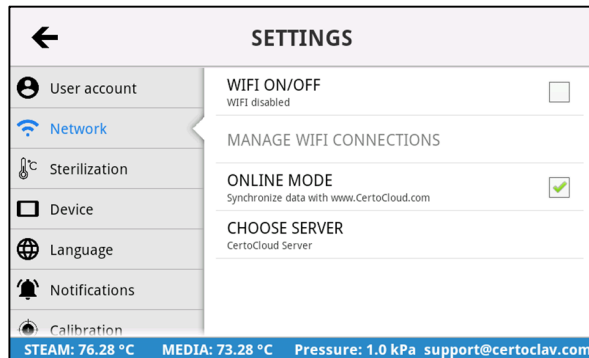

10. Select the "Add server from local network" option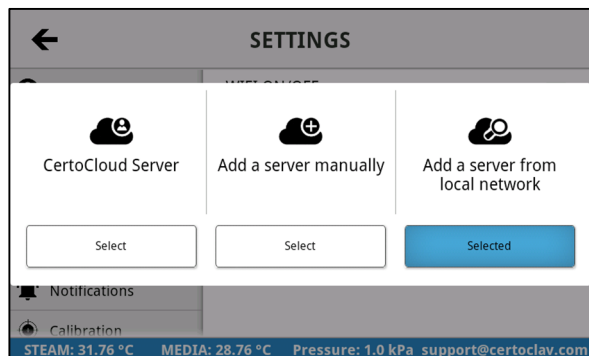


11. The autoclave searches for the CertoServer in the local network.
12. After the CertoServer has been found, confirm with the Save button.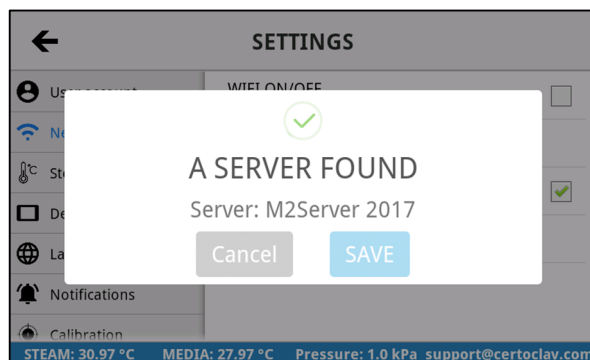

33.6 Sterilization settings

33.6.1 Export of all logs to USB stick


The touch PC can store about 1000 protocols in its internal memory. To export all protocols to USB stick, insert a USB stick into the interface box. Navigate to Settings->Sterilization->Export protocols to USB stick.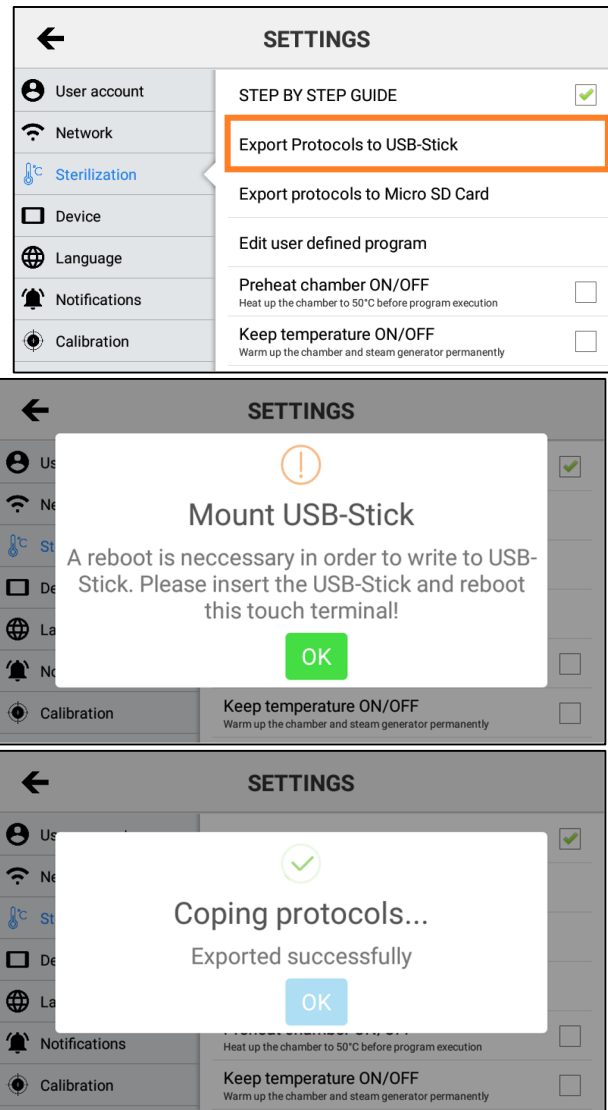

If the log export fails, your USB stick is not recognized at first attempt. In this case, leave the USB stick plugged into the device and restart the touch screen. Now try to export the logs again.


33.6.2 Export of all logs to SD card

4. Insert a micro SD card into the foldable slot on the right side of the touch screen.
5. Press the list item "Export all logs to micro SD card".
6. If the log export fails, restart the touch screen with the SD card inserted. Now try to export the logs again.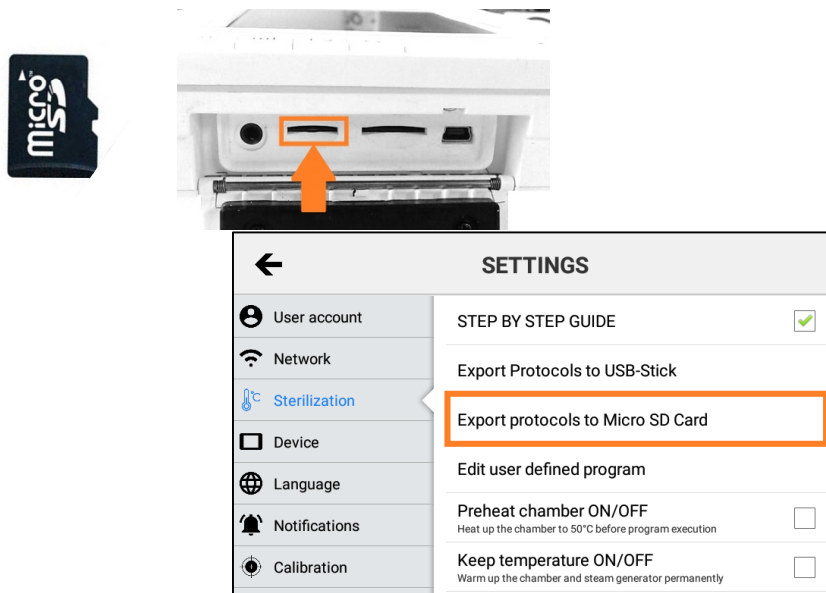


33.6.3 Export of all protocols via network

Log into the autoclave with your CertoCloud or CertoServer account. Make sure that the online mode is activated. If the login is successful, all logs are uploaded to CertoCloud. The export was successful if the Cloud icon is displayed with a check mark.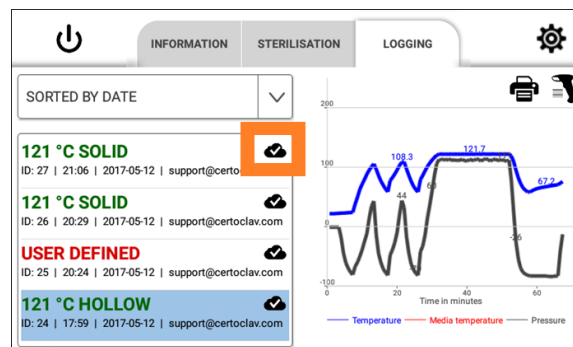

33.6.4 Label Designer


To mark items with barcodes, you can use the integrated Label Designer. The Label Designer can be opened under Settings->Sterilization->Label Designer.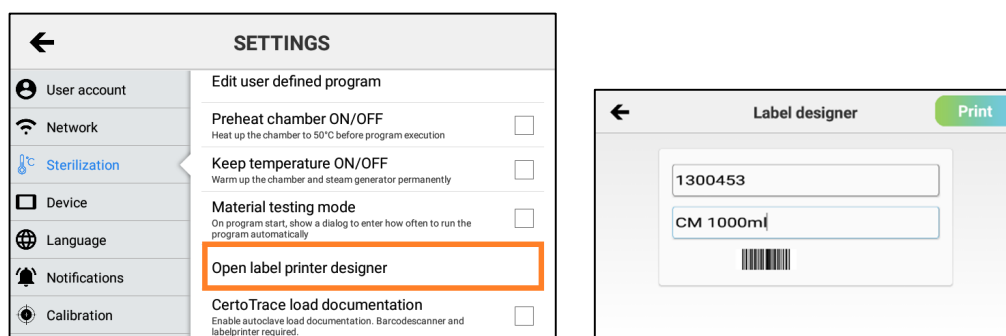

33.7 System settings


33.7.1 Install updates via Internet

5. Make sure that the autoclave is connected to the Internet.
6. Click on the "Device" entry.
7. In the details pane, scroll to the right side "Check for updates" and click on this entry.
8. A progress bar appears with the download. Once all updates have been downloaded, they will be installed automatically. If a new software version is available, an installation wizard will guide you through the installation process.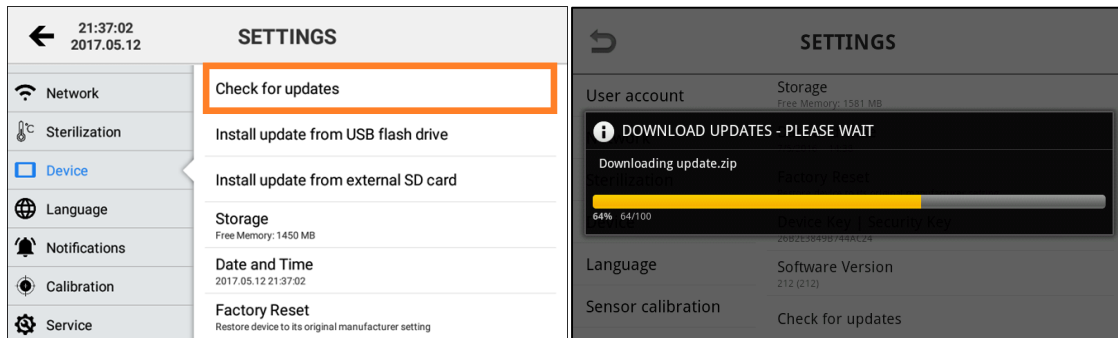


33.7.2 Install updates via USB stick

6. Download the latest update from www.certoclav.com or ask for the latest update for your device by emailing support@certoclav.com. Please provide the support team with the serial number of the autoclave. The serial number can be found on the type plate of the autoclave or under Settings->Device->Serial number. Copy the "update.zip" to the root directory of a USB stick and insert this USB stick into the interface box.
7. Click on the "Device" entry.
8. In the detail view on the right side, scroll down to the entry "Installing updates via USB stick" and press this entry.
9. If the USB stick is not recognized, restart the touch screen and try again.
10. Follow the on-screen instructions to install the appropriate updates.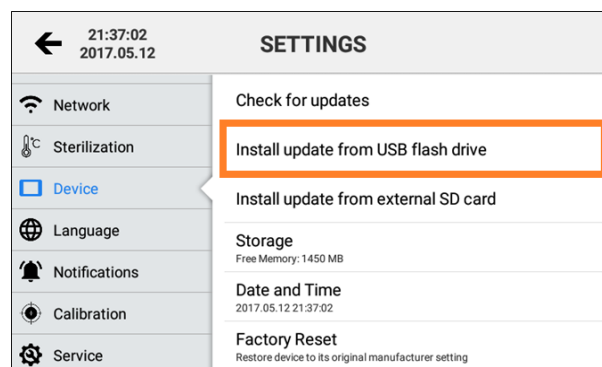

33.7.3 Install updates via SD card

- Download the latest update from www.certoclav.com or ask for the latest update for your device by e-mail to support@certoclav.com. Please provide the support team with the serial number of the autoclave. The serial number can be found on the type plate of the autoclave or under Settings->Device->Serial number. Copy the "update.zip" to the root directory of a micro SD card and insert this SD card into the slot on the right side of the touch screen.
- Click on the "Device" entry.
- In the detail view, scroll to the right side "Installing updates via SD card" and press this entry.
- If the SD card is not recognized, restart the touch screen and try again.
- Follow the on-screen instructions to install the appropriate updates.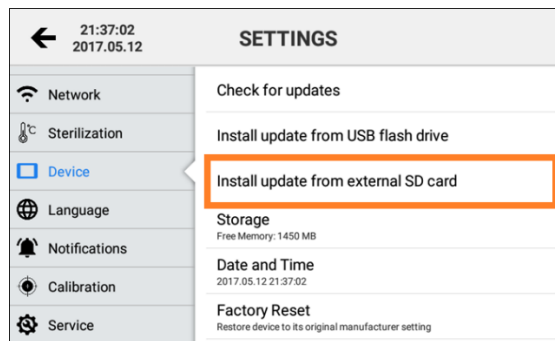

33.7.4 Restore factory settings


If you reset the touchscreen to factory settings, all logs and user accounts not previously exported and not synchronized with CertoCloud will be deleted.

- Click on the Device entry.
- Scroll to the Restore factory settings entry in the detail view on the right side and press this entry.
- The Touch PC was reset. All saved user accounts and logs have been deleted.

33.7.5 Change language

4. In the settings menu, open the language menu by clicking the language button.
5. Press current language in the detail view on the right side of the screen
6. You will see a scrollable list of supported languages. Select a language by clicking on the desired language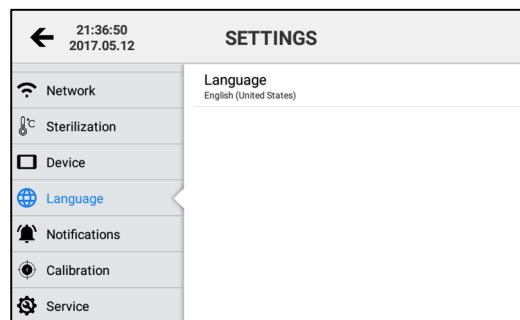


Finally, press "Save" to fix the changes.

33.7.6 Configuration of SMS and e-mail notification

If you do not have a CertoCloud account yet, you can create a new CertoCloud account at the login window. Note: To create the CertoCloud account, the autoclave must be connected to the Internet. Log in with your CertoCloud account.

Navigate  Settings -> Notifications to configure SMS and email notifications. If you want to receive an e-mail when a program has been cancelled, maintenance is required, or a program has been completed successfully, then activate the respective e-mail checkbox. Enter the email address of your choice which should receive the notification. If you have a CertoCloud Premium account, you can also enter a cell phone number that will receive a notification via SMS.

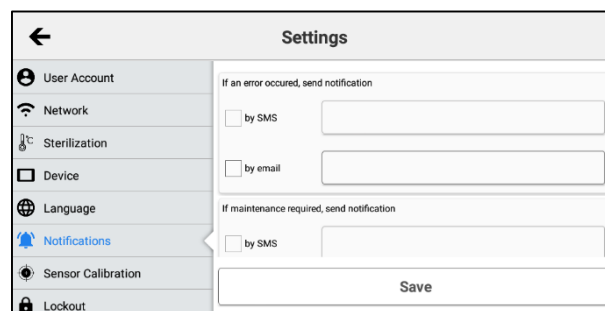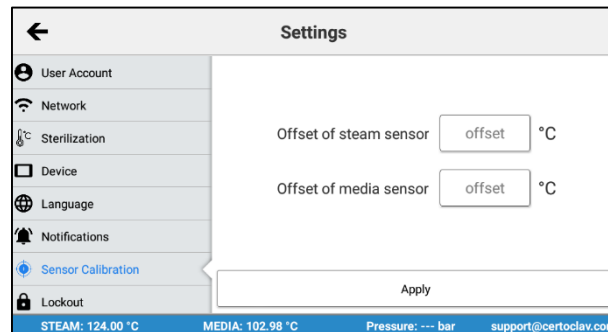
Finally, press the Save field at the bottom of the screen to save all the changes you have made.

33.8 Sensor calibration

Sensor calibration should only be performed by trained personnel.

4. Navigate to  Settings-> Calibration to display the Sensor Offset parameters.
5. Press the offset you want to adjust. The offset parameters may be in the interval [-3°C; 3°C].
6. Then press the Apply button.

33.8.1 Audit Trail

An audit trail is a quality assurance tool designed to control and record changes made to processes. With audit trails, the focus is on monitoring the change and deletion actions of the various users. Thus, interventions in processes become comprehensible, controllable and fully visible to downstream users.

Time	User	Event ID
Feb 03, 2021 02:45 PM	Admin	33
Feb 03, 2021 02:45 PM	Admin	35
Feb 03, 2021 02:44 PM	Admin	18
Feb 03, 2021 02:44 PM	Admin	18
Feb 03, 2021 02:44 PM	Admin	18
Feb 03, 2021 02:44 PM	Admin	18
Feb 03, 2021 02:44 PM	Admin	15
Feb 03, 2021 02:44 PM	Admin	15

The status bar at the bottom shows: STEAM: 89 °C, MEDIA: 86 °C, PRESS: 1 kPa, support@certoclav.com.

Subsequently, it is also possible to learn more about the event. To do this, click on a desired time and get more detailed information about the event. A dialog box will appear for this purpose.


Time	User	Event ID
Feb 03, 2021 02:44 PM	Admin	18
Old Value		New Value
preferences_list_pressure_unit = bar		preferences_list_pressure_unit = kPa

The dialog box has an 'OK' button. The background table shows the same audit logs as the previous screenshot. The status bar at the bottom shows: STEAM: 81 °C, MEDIA: 78 °C, PRESS: 1 kPa, support@certoclav.com.

33.8.2 Lock settings

To prevent unauthorized changes to the autoclaves, sensor parameters, programs and settings, the settings can only be changed in the password-protected administrator user

account. Here you can define whether other user roles have access to certain settings.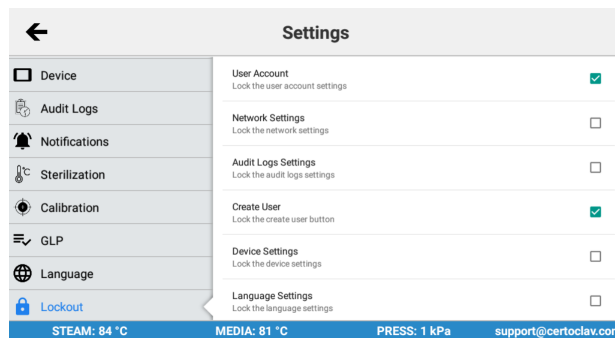


Setting	Option(default value is marked bold)
User account	YES , NO
Network settings	YES, NO
Audit logs settings	YES, NO
Create user	YES , NO
Devices settings	YES, NO
Language settings	YES, NO
Notification settings	YES, NO
Sterilization settings	YES, NO
Calibration settings	YES, NO
GLP settings	YES, NO
Service settings	YES, NO

Here "YES" means that the box is checked by default and the menu is not visible for the other user roles.

34 PC software

34.1 CertoCloud Software Features

CertoCloud is a browser-based management software for your autoclave. CertoCloud is available at www.certocloud.com for PC, smartphone and tablets. CertoCloud is also available as a lab-internal private cloud. For this, CertoServer must be connected to your laboratory network. CertoServer does not require an Internet connection. Type <http://certocloud> into your web browser to open your PC software.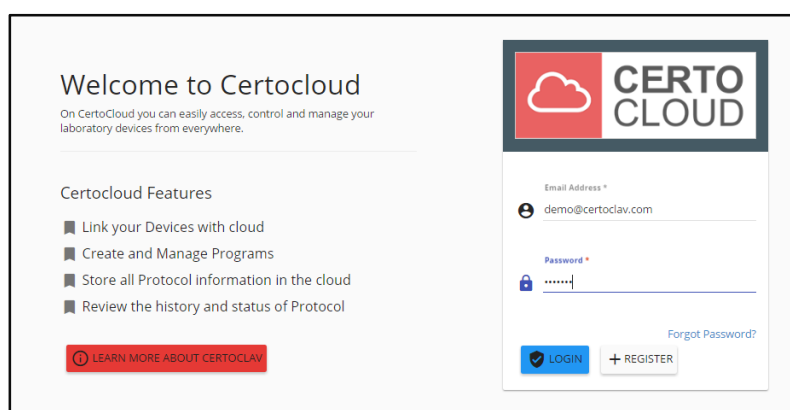

Create a user account, directly at the autoclave or web-based

Email Address *
max.muster@mail.com

Password *
.....

Forgot Password?

LOGIN + REGISTER NOW

Tracking real-time data from the autoclave

Autoclave Lab 2

Serial
220002576

Model
CertoClav
22 Vac
Pro

Firmware
DEFAULT

Management of SMS and email notifications. For example, when maintenance is due.

If An Error Occurred, Send Notification

SMS 0123456789 X


EMAIL max.muster@mail.com X

Download all sterilization protocols as signed PDF, Excel and ZIP file.

List of Protocols

Cycle ↑	Program Name	Start
76	cloud generated	7/22/ PM
75	121°C SOLID	7/22/ PM
74	Neues Programm	7/22/ PM
74	121°C SOLID	7/22/ PM

Viewing of all logs in the graphical view, as a list or detailed view.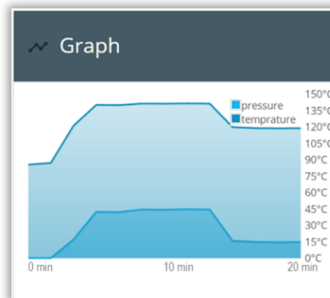

Prove the validity of your exported logs.

Protocol Verification

Protocol ID * 0 / 24


Device Key * 0 / 17

Security Code * 0 / 40

VERIFY

34.2 Display of sterilization protocols via web browser

If your autoclave is connected to the Internet or to the local CertoServer. Using a browser on the same network, visit www.certocloud.com or <http://certocloud> to manage the autoclave and view logs. An example log is shown below.

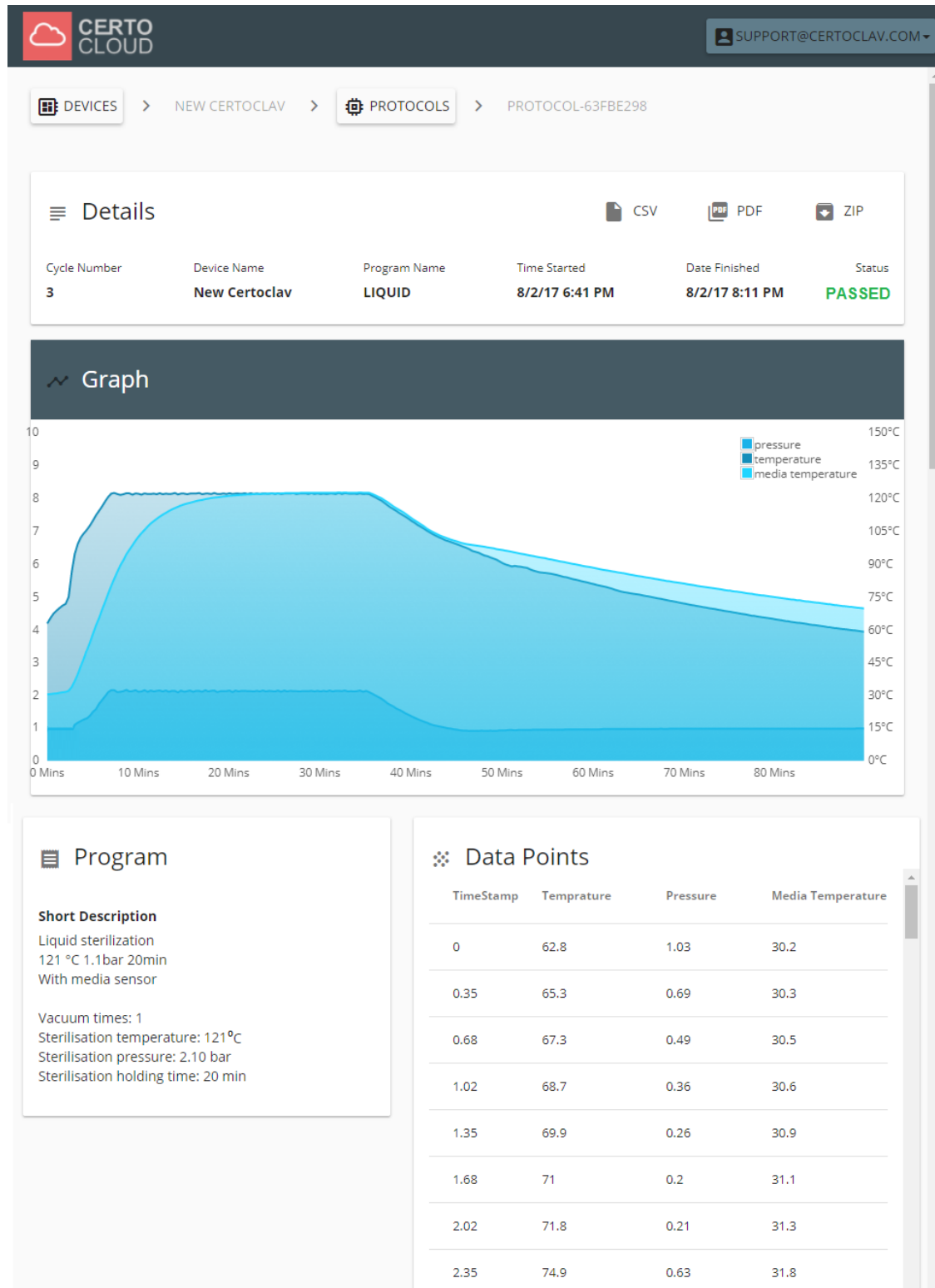35 Transport and storage

35.1 Storage conditions

The CertoClav Connect should be stored at a room temperature between 5°C and a maximum of 40°C. The relative humidity must not be higher than 80%. Ensure that the storage facility is well ventilated and does not contain any gases that promote corrosion.

35.2 Transport conditions

Make absolutely sure that all water has been drained from the sterilization chamber and the condensate container before transport.

Do not drop the CertoClav Connect during transport.

36 Warranty

Warranty: 2 years. The right to free service expires if one of the following conditions situations has occurred:

- The damage was caused by incorrect installation and/or operation.
- The damage was caused by negligent behavior.
- The damage was caused by force majeure such as unacceptable supply voltage or fire.